INCLUDED PARTS:

Before the assembly, please make sure that all the following parts are included in the package:

- Backrest, metal plate, small knob - 1 set
- Seat with metal frame - 1 set
- Long mechanism with one button - 1 set
- Column - 1 piece
- Extension of the column - 1 piece
- Footrest - 1 piece
- Column cover - short (A) and long (B) - 2 pieces
- A base of five branches - 1 piece
- Landslides - 5 pieces

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1. Fix the sliders (1) on the base of five branches
2. Fix the extension of the column (3A) to the hole in the base (2).
3. Place the column cover on the extension of the column (3A).
4. Attach the footrest (4) to the extension of the column (3A).

When mounting a footrest, the chrome edge should be placed with the black 3-arm support (on top).

5. Fix the column (3B) in the extension hole of the column (3A).
6. Insert the seat back fastening element (6) into the guides on the seat leaf (5) and secure it with a longer knob.
7. Unscrew the smallest knob on the backrest (7), insert the seat back fastening element (6) and screw the knob one more time.
8. Backrest (7) mounted with a seat (5) placed on the pin of the column (3B).
9. Adjust the height of the seat and the backrest and fix the adjustment screws.

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