



INCLUDED PARTS:

Before the assembly, please make sure that all the following parts are included in the package:

- Seat with metal plate - 1 piece
- Column - 1 piece
- Footrest with knob - 1 piece
- A base of five branches - 1 piece
- Landslides- 5 pieces

ASSEMBLY INSTRUCTIONS

1. Attach the sliders (1) to the base of five branches (2).
ATTENTION! The landslides must already be fixed at the base of five branches.

2. Fix the column (3) in the hole in the base (2).

3. Place the footrest (4) on the column (3).

ATTENTION!

When mounting a footrest, the chrome edge should be placed with the black 3-arm support (on top).

4. Place the seat (5) on the column pin (3).

5. Adjust the height of the stool.





LISTA DE PARTES INCLUIDAS:

Antes del montaje, asegúrese de que todas las piezas especificadas a continuación estén incluidas en un paquete:

- Asiento con placa de metal – 1 pieza
- Columna - 1 pieza
- Reposapiés con pomo - 1 pieza
- Una base de cinco ramas - 1 pieza
- Deslizamientos- 5 piezas

INSTRUCCIONES DE MONTAJE

1. Fije los deslizadores (1) en la base de cinco ramas (2).
¡ATENCIÓN! Los deslizamientos deben estar ya fijados en la base de cinco ramas.

2. Fije la columna (3) en el orificio de la base (2).

3. Coloque el reposapiés (4) en la columna (3).

¡ATENCIÓN!

Montando un reposapiés, el borde cromado debe colocarse en un soporte negro de 3 brazos (en la parte superior).

4. Coloque el asiento (5) en el pasador de la columna (3).

5. Ajuste la altura del taburete.

