## MM1500 Microplate Mixer

*Please read the User Manual carefully before use, and follow all operating and safety instructions!*

# user manual

english / español / français

# User manual

## MM1500 Microplate Mixer

### Preface

Users should read this Manual carefully, follow the instructions and procedures, and beware of all the cautions when using this instrument.

### Service

If help is needed, you can always contact your dealer or Labbox via [www.labbox.com](http://www.labbox.com)

Please, provide the customer service representative with the following information:

- Serial number (on the back side)
- Description of the problem
- Your contact information

### Warranty

This instrument is guaranteed to be free from defects in materials and workmanship under normal use and service, for a period of 24 months from the date of invoice. The warranty is extended only to the original purchaser. It shall not apply to any product or parts which have been damaged on account of improper installation, improper connections, misuse, accident or abnormal conditions of operation.

For claim under the warranty, please contact your supplier.

# 1. Safety Instructions

	<b>Warning!</b> <ul style="list-style-type: none"><li>• Read the operating instructions carefully before use.</li><li>• Ensure that only trained staff works with the instrument.</li></ul>
	<b>Protective ground contact!</b> <ul style="list-style-type: none"><li>• Make sure that socket must be grounded (protective ground contact) before use.</li></ul>

Table 1

- When working, wear the necessary personal protective equipment to avoid the risk of:
  - Burns caused by splashing and evaporation of liquids
  - Intoxication caused by release of toxic or flammable gases.
- Set up the instrument on a spacious, stable, clean, non-slip, dry, and fireproof surface that can support the weight of the equipment. Do not operate the instrument in explosive atmospheres, with hazardous substances, or under water.
- Gradually increase the speed. Reduce the speed if:
  - The medium splashes out of the vessel because of too high speed.
  - The appliance is not running smoothly.
- Beware of hazards due to:
  - Overfilling of media
  - Unsafe container
- Check the appliance and accessories beforehand for damage each time you use them. Do not use damaged components. Safe operation is only guaranteed with the accessories described in the “Accessories” chapter. Accessories must be securely attached to the device and cannot come off by themselves.
- Always disconnect the plug before assembling or disassembling accessories.
- The equipment can only be disconnected from the power supply by pulling out the mains power cable.
- Make sure the power supply voltage match the power needs of the equipment.
- The instrument may only be opened by experts.
- Keep away from high magnetic fields.

- Observe the minimum distances between devices, between the device and the wall, and above the assembly (min. 100 mm).

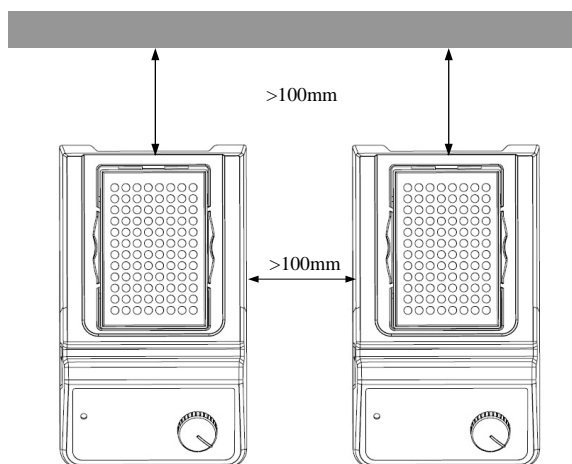
Figure 1

## 2. Proper Use

The instrument is designed for mixing liquids in schools, laboratories, or factories. This device is not suitable for use in residential areas or other restrictions mentioned in Chapter 1.

## 3. Inspection

### 3.1 Unpacking

Unpack the equipment carefully and check for any damages that may have arisen during transportation. Please contact the manufacturer/supplier for technical support.

**Note:**

If there is any apparent damage to the system, please do not plug it into the power line.

### 3.2 Listing of Items

The package includes the following items:

Items	Qty
Main unit (with adapter for 1 microplate)	1
Power cable	1
User Manual	1

Table 2

## 4. Control

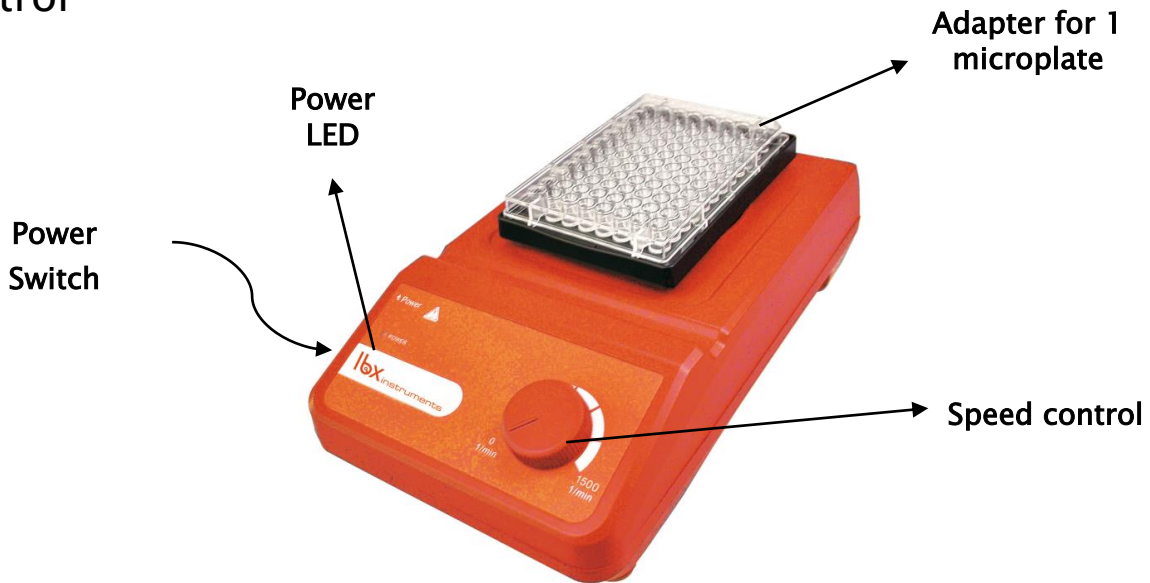
Figure 2 : MM1500 series

	Items	Descriptions
MM1500	Speed control knob	The function “mixing” is switched ON or OFF with the knob. To increase the value, turn the knob clockwise.
	Power Switch I/O	Switch the instrument ON or OFF.
	Power LED	While the device is ON, the power LED is lit.

Table 3

## 5. Trial Run

- Make sure that the required operating voltage and power supply voltage match.
- Ensure the socket is properly earthed.
- Plug in the power adapter; ensure the power is on.
- Switch ON the instrument by pressing the power switch I/O.
- Set the mixing speed with the control knob. The device begins to work.
- Switch OFF the instrument by pressing the power switch I/O.

*If these operations above cause no abnormalities, the device is ready to operate. Otherwise, the device may have been damaged during transportation; please contact your manufacturer/supplier for technical support.*

### Warning!

Forbidden to transfer the vessel when the instrument is working.

## 6. Operation

### 6.1 Placing and removing the microplate

The microplate is to be placed properly, holding it from one of the sides as shown in figure 3.

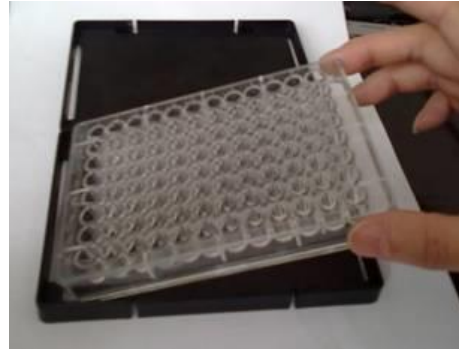
Figure 3

### 6.2 Speed range

- Microplate clamp: 0–1500rpm.
- Double microplate clamp: 0–1000rpm.

## 7. Maintenance and Cleaning

- Proper maintenance can keep instruments working properly and lengthen their lifetime.
- Do not spray cleanser onto the instrument when cleaning.
- Unplug the power line when cleaning.
- Only use recommended cleansers:

Dyes	Isopropyl alcohol
Construction materials	Water containing surfactants / Isopropyl alcohol
Cosmetics	Water containing surfactants / Isopropyl alcohol
Foodstuffs	Water containing surfactants
Fuels	Water containing surfactants

Table 4

- Before using any other cleaning or decontamination methods, verify with the manufacturer that such method will not harm the equipment.
- Keep the device clean and protect it from liquid splashes, which could shorten its lifetime.
- Wear the proper protective gloves while cleaning the instrument.

**Note:**

- Electronic devices cannot be cleaned with cleanser.
- If maintenance service is needed, the device must be cleaned before shipping to avoid pollution from hazardous substances, and shipped in its original packaging.
- If the instrument will not be used for a long time, please switch it off and place it in a dry, clean, and stable location at room temperature.

## 8. Standards and Regulations

Manufacture in accordance with the following safety standards:

EN 61010-1

Manufacture in accordance with the following EMC standards:

EN 61326-1

Associated EU guidelines:

EMC-guidelines: 89/336/EWG

Machine guidelines: 73/023/EWG

## 9. Specifications

Items	Specifications
Voltage [VAC]	100~240
Frequency [Hz]	50/60
Power [W]	20
Movement type	Orbital
Orbital diameter [mm]	4.5
Max. mixing weight (with attachments)	0.5
Motor type	Exterior rotor Brushless motor
Motor rating input [W]	18
Motor rating output [W]	10
Permissible ON time [%]	100
Speed range [rpm]	0 - 3000
Speed display	Scale
Microplate working	Yes
Microplate amount	1
Dimensions (mm)	260×150×80
Weight [kg]	3
Permitted ambient temperature[°C]	5-40
Permitted relative humidity	80%
Protection class acc. to DIN 60529	IP21

Table 5

## 10. Accessories

Please contact our company to order the following accessories available for this device:

Accessories	References	Descriptions	Running mode
	MIXS-A31-001	Universal top plate (round). Used with tube adapters	Continuous operation
	MIXS-A11-001	Adapter for 1 microplate	Continuous operation
	MIXS-A21-001	Adapter for 2 microplates	Continuous operation
	MIXS-A41-001	Tube adapter for 48 hole test tubes, Ø 6 mm, used with MIXS-A31-001	Continuous operation
	MIXS-A51-001	Tube adapter for 18 hole test tubes, Ø 10 mm, used with MIXS-A31-001	Continuous operation
	MIXS-A61-001	Tube adapter for 12 hole test tubes, Ø 12 mm, used with MIXS-A31-001	Continuous operation
	MIXS-A71-001	Tube adapter for 8 hole test tubes, Ø 16 mm, used with MIXS-A31-001	Continuous operation
	MIXS-A81-001	Tube adapter for 8 hole test tubes, Ø 20 mm, used with MIXS-A31-001	Continuous operation

Table 6

Changing accessories:

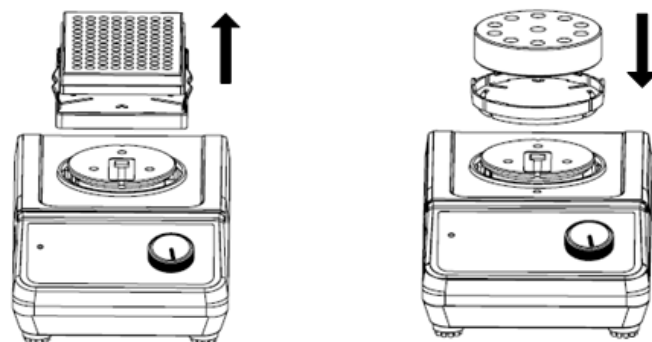
Figure 4

**Remarque:**

Be careful when replacing accessories, otherwise over time these accessories may be damaged.

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